

RT-qPCR Lyo-Ready Mix FOR TRIAL

Description

RT-qPCR Lyo-Ready Mix FOR TRIAL is a universal one-step probe product designed for evaluating the RT-qPCR Lyo-Ready Mix (#4501) performance before committing to large-scale lyophilization projects. This trial version is intended for performance testing only and is not fully lyophilization-ready.

Unlike the standard RT-qPCR Lyo-Ready Mix (#4501), which includes the 10,000× lyophilization-ready reverse transcriptase (RTase Lyo) and two dilution buffers, this trial kit provides the RTase Lyo in a diluted 20x format, well suited for performance testing. It delivers the same robust, sensitive, and fast RT-qPCR performance, but is not suitable for lyophilization.

If lyophilization studies are required, we recommend using the RT-qPCR Lyo-Ready Mix (#4501), which is specifically optimized for that application.

The mix features state-of-the-art technology, including an antibody-regulated hot-start Taq polymerase, optimized buffer chemistry, PCR enhancers, RNase inhibitor and stabilizers, enabling efficient cDNA synthesis and real-time PCR amplification in a single tube.

Compatible with a range of probe chemistries such as TaqMan® and Scorpions®, this trial mix supports rapid detection and quantification of diverse RNA targets, including mRNA, viral RNA, and total RNA.

This trial format is supplied in a size suitable for evaluation (125 reactions) making it easier and more cost-effective to assess assay performance before scaling up for lyophilization testing.

Kit Components

Component	Volume
RT-qPCR Lyo-Ready Mix (2x)	1 x 1.25 mL
RTase Lyo (20x)	1 x 0.125 mL

Storage and Shipment

Transport with an ice pack. The reagents should be stored at -20°C upon arrival. The reagents are stable until the expiration date if stored correctly.

Reaction Master Mix Set-Up

The recommended master mix set-up for a 20 µL reaction volume is shown in the table below.

Reagent	Volume (µL)	Final concentration
RT-qPCR Lyo-Ready Mix (2x)	10	1x
∞Forward primer (10 µM)	X	100–500 nM
∞Reverse primer (10 µM)	X	100–500 nM
∞Probe (10 µM)	X	50–250 nM
RTase Lyo (20x)	1	1x
RNA template	2–8	Variable
Nuclease-free Water	Up to 20 µL final volume	

∞Primers and probes should be specific to the target DNA/RNA of interest. The recommended T_m for primers is between 56°C and 60°C, and the T_m for probes should be between 65°C and 70°C.

Instrument and Program Set-Up

Cycles	Steps	Temperature	Time
1	Reverse transcription	50°C	10 min
1	Polymerase activation	95°C	1 min
40	Denaturation	95°C	5 sec
	Annealing/Extension	60°C	20 sec



Legal disclaimer

Technical information and support

For technical enquiries or assay development support, please contact us via e-mail at: mdx@medixbiochemica.com. Additional information and technical resources are available on our website at: info.medixbiochemica.com/resources.



Legal disclaimer

myPOLS Biotec part of Medix Biochemica group products meet their specifications if transported, stored, and used according to the instructions. Medix Biochemica's products may not be used or reproduced without Medix Biochemica's written permission.